ACCELERATED SURFACE AREA AND POROSIMETRY SYSTEM



micromeritics[®]

Effective Solutions for Material Characterization

CALCULATIONS



CORPORATE PROFILE

Micromeritics Instrument Corporation is a leading global provider of solutions for material characterization with best-in-class instrumentation and application expertise in five core areas: density; surface area and porosity; particle size and shape; powder characterization; and catalyst characterization and process development. Founded in 1962, the company is headquartered in Norcross, Georgia, USA and has more than 300 employees worldwide. With a fully integrated operation that extends from a world class scientific knowledge base through to in-house manufacture, Micromeritics delivers an extensive range of high-performance products for academic research and industrial problem-solving. The implementation of tactical partnerships to incubate and deliver valuable new technologies exemplifies the company's holistic, customer-centric approach which extends to a cost-efficient contract testing laboratory – the Particle Testing Authority (PTA). The strategic acquisitions of Freeman Technology Ltd and Process Integral Development S.L. (PID Eng & Tech) reflect an ongoing commitment to optimized, integrated solutions in the industrially vital areas of powders and catalysis.

Freeman Technology (Tewkesbury, UK) brings market-leading powder characterization technology to Micromeritics' existing portfolio of particle characterization techniques. The result is a suite of products that directly supports efforts to understand and engineer particle properties to meet powder performance targets. With over 15 years of experience in powder testing, Freeman Technology specializes in systems for measuring the flow properties of powders. In combination with detailed application know-how these systems deliver unrivalled insight into powder behavior supporting development, formulation, scale-up, processing and manufacture across a wide range of industrial sectors.

PID Eng & Tech (Madrid, Spain) complements Micromeritics' renowned offering for catalyst characterization with technology for the measurement and optimization of catalytic activity, with a product range that extends to both standard and bespoke pilot scale equipment. Launched in 2003, PID Eng & Tech is a leading provider of automated, modular microreactor systems for the detailed investigation of reaction kinetics and yield. These products are supported by a highly skilled multidisciplinary team of engineers with in-depth expertise in the design, construction and operation of laboratory units and process scale-up.

The Particle Testing Authority (PTA) provides material characterization services for fine powders and solid materials using Micromeritics' instrumentation alongside complementary solutions from other vendors. With the certification and expertise to operate across a wide range of industries the PTA offering runs from single sample analysis to complex method development, method validation, new product assessment, and the analytical support required for large scale manufacturing projects. An experienced, highly trained team of scientists, engineers, and lab technicians works closely with every client to ensure that all analytical requirements are rapidly and responsively addressed.

Micromeritics has a strong global network with offices across the Americas, Asia, and Europe complemented by a dedicated team of distributors in additional locations. This ensures that local, knowledgeable support is available for every customer, in academia or industry. Micromeritics works across a truly diverse range of industries from oil processing, petrochemicals and catalysts, to food and pharmaceuticals, and at the forefront of characterization technology for next generation materials such as graphene, metal-organic-frameworks, nanocatalysts, and zeolites. Engineering solutions that work optimally for every user is a defining characteristic of the company.



CONTACT US

Micromeritics Instrument Corporation

4356 Communications Drive Norcross, GA / USA / 30093-2901

Phone: 1-770-662-3636 Fax: 1-770-662-3696 www.Micromeritics.com

Instrument Service or Repair

Phone: 1-770-662-3666

International — contact your local distributor or call 1-770-662-3666

Service.Helpdesk@Micromeritics.com

Micromeritics Learning Center

Phone: 1-770-662-3607

www.Micro.edu



Table of Contents

Corporate Profile	ii
Contact Us	iii
Alpha-S Method	1
BET Surface Area	2
BJH Pore Volume and Area Distribution	3
Crystallite Size for Chemical Adsorption	12
Data Smoothing for Chemical Adsorption	13
DFT (Density Functional Theory)	14
Dollimore-Heal Pore Volume and Area Distribution	17
Dubinin-Astakhov	19
Dubinin-Radushkevich	23
Equation of State	25
Equilibration	26
f-Ratio Method	27
Free Space for Chemical Adsorption	28
Freundlich Isotherm	29
Heat of Adsorption for Physical Adsorption	30



lorvath-Kawazoe	31
Langmuir Surface Area for Chemical Adsorption	41
Langmuir Surface Area for Physical Adsorption	42
Metal Dispersion for Chemical Adsorption	44
Metallic Surface Area	45
MP-Method	46
Quantity Adsorbed for Chemical Adsorption	48
Quantity Adsorbed for Chemical Adsorption	49
Quantity Adsorbed for Physical Adsorption	50
Relative Pressure	52
Saturation Pressure	53
SPC Report Variables	54
Stoichiometry Factor	56
Summary Report	57
t-Plot	59
Temkin Isotherm	60
Thermal Transpiration Correction	61
Thickness Curve	62



ALPHA-S METHOD

The alpha-S curve is calculated from the reference isotherm by dividing each quantity adsorbed by the quantity adsorbed at 0.4 relative pressure.

$$lpha_i = rac{Q_i}{Q_{0.4}}$$

where $Q_{0.4}$ is found by linear interpolation.

A least-squares analysis fit is performed on the $(\alpha_i, Q_{ads,i})$ data pairs where α_i is the independent variable and $Q_{ads,i}$ is the dependent variable. The following are calculated:

- a. Slope (S cm³/g STP)
- b. Y-intercept (Q₀ cm³/g STP)
- c. Error of the slope (cm³/g STP)
- d. Error of the Y-intercept (cm³/g STP)
- e. Correlation coefficient

Surface area is calculated as:

$$A_s=rac{A_{
m ref}S}{Q_{0.4}}$$

Pore size is calculated as:

$$rac{Q_0V_{
m mol}}{22414{
m cm}^3{
m STP}}$$

where $V_{
m mol}$ is liquid molar volume from the fluid property information.



BET SURFACE AREA

The BET¹) transformation is calculated as:

$$y=rac{1}{Q(
ho^0/
ho-1)}$$

A least-squares fit is performed on the $P_{
m rel}$, ${f y}$. The following are calculated:

- a. Slope (S cm³/g STP)
- b. Y-intercept $(y_0 \text{ cm}^3/\text{g STP}))$
- c. Uncertainty of the slope $(u(s) cm^3/g STP))$
- d. Uncertainty of the Y-intercept $(u(y_0) cm^3/g STP)$
- e. Correlation coefficient

Using the results of the above calculations, the following can be calculated:

BET Surface Area (m²/g):

$$A_s = rac{A_m N_A}{V_m (s + y_0)} imes 10^{-18} m^2 / nm^2$$

BET C value:

$$C = rac{s}{y_0} + 1$$

Quantity in the Monolayer (cm³/g STP):

$$Q_m=rac{1}{Cy_0}=rac{1}{s+y_0}$$

Error of the BET Surface Area (m²/g):

$$u(A_s)=rac{\sqrt{u^2(s)+u^2(y_0)}}{s+y_0}$$

¹⁾ Brunauer, S.; Emmett, P.H.; and Teller, E., J. Am. Chem. Soc. 60, 309 (1938).



BJH PORE VOLUME AND AREA DISTRIBUTION

For adsorption data, the relative pressure and quantity adsorbed data point pairs collected during an analysis must be arranged in reverse order from which the points were collected during analysis. All calculations are performed based on a desorption model, regardless of whether adsorption or desorption data are being used.

The data used in these calculations must be in order of strictly decreasing numerical value. Points which do not meet this criterion are omitted. The remaining data set is composed of relative pressure (P), quantity adsorbed (Q) pairs from (P_1, Q_1) to (P, Q_n) where $(P_n = 0, Q_n = 0)$ is assumed as a final point. Each data pair represents an interval boundary (or desorption step boundary) for intervals i=1 to i=n-1 where n = total number of (P, Q) pairs.

Generally, the desorption branch of an isotherm is used to relate the amount of adsorbate lost in a desorption step to the average size of pores emptied in the step. A pore loses its condensed liquid adsorbate, known as the core of the pore, at a particular relative pressure related to the core radius by the Kelvin¹) equation. After the core has evaporated, a layer of adsorbate remains on the wall of the pore. The thickness of this layer is calculated for a particular relative pressure from the thickness equation. This layer becomes thinner with successive decreases in pressure, so that the measured quantity of gas desorbed in a step is composed of a quantity equivalent to the liquid cores evaporated in that step plus the quantity desorbed from the pore walls of pores whose cores have been evaporated in that and previous steps. Barrett, Joyner, and Halenda²) developed the method (known as the BJH method) which incorporates these ideas. The algorithm used is an implementation of the BJH method.

EXPLANATION OF TERMS


A pore filled with condensed liquid nitrogen has three zones:

- Core. The core evaporates all at once when the critical pressure for that radius is reached; the relationship between the core radius and the critical pressure is defined by the Kelvin equation.
- Adsorbed layer. The adsorbed layer composed of adsorbed gas that is stripped off a bit at a time with each pressure step; the relationship between the thickness of the layer and the relative pressure is defined by the thickness equation.
- Walls of the cylindrical pore. The walls of the cylindrical pore the diameter of the empty pore is required to determine the pore volume and pore area. End area is neglected.

¹⁾ Kelvin, J. (published under the name of Sir William Thomson), Phil. Mag. 42, 448-452 (1871).

²⁾ Barrett, E.P.; Joyner, L.S.; and Halenda, P.P., J. Am. Chem. Soc. 73, 373-380 (1951).



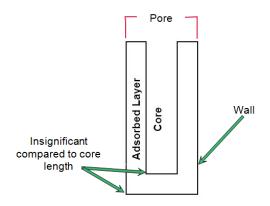

CALCULATIONS

The quantities adsorbed (Qa) are converted to the liquid equivalent volumes (VI, cm³/g)

$$V l_i = rac{Q_i V_{
m mol}}{22414 {
m cm}^3 {
m STP}}$$

where $V_{\rm mol}$ is the liquid molar volume from the fluid property information.

The relative pressure (P_i) is assumed to be close to unity so that substantially all the pores in the sample are filled.

The corresponding Kelvin core radius is calculated. Only pores smaller than this size will be included:

$$Rc_i = rac{-A}{(l+F)\ln(P_i)}$$

where

A = adsorbate property factor (from the BJH Adsorptive Options window)

F = fraction of pores open at both ends (from the BJH Adsorption Report Options window or the BJH Desorption Report Options window); assumed to be zero for desorption

Rc = Kelvin radius (Å) of core



This radius will be adjusted for the thickness of the adsorbed layer during subsequent calculation steps

The following calculations (a-c) are made for each relative pressure interval based on the increment of volume desorbed during that interval:

where

- i = interval number, that is i=1 for the first interval from P1 to P2, and so on
- j = each previous interval during which new pores were found
- k = the total number of intervals in which new pores have been found. It is also the number of lines reported on the BJH table for collected data
- a. The thickness of the adsorbed layer at the end of the interval is calculated using the equation located in *Thickness Curve on page 62*.

For the last pressure interval from the lowest Pr_i to zero relative pressure, reference the calculations from the equations in *Thickness Curve on page 62*.

For the first pressure interval, there are no previously opened pores so the volume of liquid desorbed from walls of previously opened pores is zero ($Vd_1 = 0$), and the remainder of Step (a) is skipped.

The change in thickness of the wall layer due to desorption from previously opened pores is calculated as:

$$\Delta Tw = Tw_1 - Tw_{i+1}$$

The annular cross-sectional area of the wall layer desorbed is calculated for all previously opened pores:

$$CSA_{
m j} = \piigg[igg(Rc_{
m j} = \Delta Twigg)^2 - Rc_{
m j}^2igg]igg(10^{-16}rac{
m cm^2}{{
m \dot a}^2}igg)$$

The total volume of gas desorbed from walls of previously opened pores is calculated:

$$Vd_{f i} = rac{\sum\limits_{f i} (LP_{f j})(CSAa_{f j})}{}$$
 for all previously opened pores

where LP_{j} = length of previously opened pores as calculated in Step b(2).



- b. The physical processes occurring for this pressure interval are determined as:
 - 1. If Vd_i is greater than the current increment of volume desorbed $(Vl_i Vl_{i+1})_{,}$ desorption from walls only is occurring. Total surface of walls exposed thus far (cm²/g) is calculated as:

$$SA_{
m w} = rac{\sum\limits_{
m j} \pi(LP_{
m j})(D_{
m avg,j})(rac{10^{-8}{
m cm}}{
m \dot{A}})}{
m eta}$$
 for all previously opened pores

where

 $D_{\text{avg.i}}$ = weighted average pore diameter calculated in Step b.2.

A new layer thickness (ΔTw) that will not overcompensate for the actual volume desorbed in this interval is calculated:

$$\Delta Tw = rac{(Vl_{ ext{i}} - Vl_{ ext{i+1}})(10^8 rac{\dot{ ext{A}}}{ ext{cm}})}{SAw_{ ext{i}}}$$

Since no cores are evaporated in this pressure interval, no new pores are revealed. Thus no ending Kelvin radius and average pore diameter are calculated for this interval. Note that this means the report may have fewer tabulated intervals on the collected data report than experimental pressure intervals.

2. If Vd_i is less than the volume increment desorbed during this interval $(Vl_i - Vl_{i+1})$, the remaining volume is due to new pores with core evaporation taking place in this interval. K, the number of intervals with new pores exposed, is increased by 1. (For the interval from the lowest Pr_1 to zero relative pressure, no new pore volume is calculated and the rest of Step b is skipped.)

The volume desorbed from newly opened pores in this interval is calculated as:

$$Vc_{\mathrm{i}} = (Vl_{\mathrm{i}} - Vl_{\mathrm{i+1}}) - Vd_{\mathrm{i}}$$

The Kelvin radius for the end of the interval is calculated as:

$$Rc_{k+1} = \frac{-A}{(1+F)\ln(P_{i+1})}$$

All new pores opened in this interval are represented by one pore having a length-weighted average pore diameter and a corresponding length sufficient to account for the required volume of adsorbate. The weighted average pore diameter is calculated as:

$$D_{
m avg,k} = rac{2(Rc_k + Rc_{k+1})(Rc_k)(Rc_{k+1})}{Rc_k^{2} + Rc_{k+1}^{2}}$$

 $D_{avg,k}$ is the diameter of a pore which would have a surface area that is the average of the areas for pores radius Rc_k and Rc_{k+1} , if its length was the mean of the lengths at those radii.



The relative pressure corresponding to D_{avg,k} is calculated as:

$$P_{ ext{avg,k}} = \expigg[rac{-2A}{(1+F)(D_{ ext{avg,k}})}igg]$$

The thickness of the adsorbed layer at this pressure is calculated as:

$$Tw_{ ext{avg,k}} = HP1igg[rac{HP2}{ln(P_{ ext{avg,k}})}igg]^{HP3}$$

The decrease in thickness of the wall layer by desorption from the walls of new pores during the lower portion of the pressure interval is calculated as:

$$\Delta Td = Tw_{ ext{avg,k}} - Tw_{ ext{i+1}}$$

The cross-sectional area of the newly opened pores is calculated as:

$$CSAc_{ extsf{k}} = \left[rac{D_{ ext{avg,k}}}{2} + \Delta Td
ight]^2 \left(rac{10^{-16} ext{cm}^2}{ ext{Å}^2}
ight)$$

The length of the newly opened pores is calculated as:

$$LP_{\mathbf{k}} = \frac{Vc_{\mathbf{i}}}{CSAc_{\mathbf{k}}}$$

Pore diameters and radii are adjusted for the change in thickness of the adsorbed wall layer during this interval. If new pores were opened during this interval, the average diameter is adjusted by the change in layer thickness during the second portion of the desorption interval as:

$$D_{ ext{avg,k,new}} = D_{ ext{avg,k,old}} + 2(\Delta T d)$$

The layer thickness change during the whole interval is added to diameters of previously opened pores as:

$$D_{ ext{avg,k,new}} = D_{ ext{avg,k,old}} + 2(\Delta T dw)$$

(not including Davg,k)



The layer thickness change desorbed during this interval also is added to the radii corresponding to the ends of the pressure intervals as:

$$Rc_{ ext{j,new}} = Rc_{ ext{j,old}} + \Delta Tw$$

for all except Rck+1.

Steps a to c are repeated for each pressure interval.

After the above calculations have been performed, the diameters corresponding to the ends of the intervals are calculated as:

$$Dp_{
m j}=2(Rc_{
m j})$$

for all Rc_i. including Rc_{k+1}.

The remaining calculations are based on Dp_i , $D_{avg,i}$, and LP_i . These calculations are only done for $D_{avg,i}$ values that fall between the Minimum BJH diameter and the Maximum BJH diameter specified by the operator on the *BJH Adsorption Report Options* window or the *BJH Desorption Report Options* window.

(1) Incremental Pore Volume (Vp_i, cm³/g):

$$Vp_{
m i}=\pi(Lp_{
m i})iggl[rac{D_{
m avg,i}}{2}iggr]^2iggl[rac{10^{16}{
m cm}^2}{{
m \dot{A}}^2}iggr]$$

(2) Cumulative Pore Volume (VP_{cum.i}, cm³/g):

$$VP_{ ext{cum,i}} = rac{\sum}{j} VP_{ ext{j}} \; ext{ for } \; (J \leq 1)$$

(3) Incremental Surface Area SA_{cum,i},m²g-A):

$$SA_{
m i} = \piigg(LP_{
m i}igg)igg(rac{10^{-2}{
m m}}{
m cm}igg)igg(D_{
m avg,i}igg)igg(rac{10^{-10}{
m m}}{
m \dot{A}}igg)$$

(4) Cumulative Surface Area SA_{cum,i},m²g):

$$SA_{ ext{cum},10} = \sum SA_{ ext{j}} ext{for} J \leq 1$$

(5) dV/dD pore volume $(dV/dD_i, cm^3/g-A)$:

$$rac{dV}{dD_{
m i}} = rac{VP_{
m i}}{Dp_{i}-Dp_{
m i+1}}$$



(6) $dV/d \log (D)$ pore volume $(dV/d \log(D)_i, cm^3g)$:

$$rac{dDv}{d\log D_{
m i}} = rac{VP_{
m i}}{\log\left(rac{Dp_{
m i}}{Dp_{
m i+1}}
ight)}$$

(7) dA/dD pore area $(dA/dD_i, m^2/g-A)$:

$$rac{dA}{dD\mathrm{i}} = rac{SA_\mathrm{i}}{Dp_\mathrm{i} - Dp_\mathrm{i+1}}$$

(8) $dA/d \log(D)$ pore area $[dA/d \log(D)i, m^2/g]$:

$$rac{dA}{d\log D_{
m i}} = rac{SA_i}{\log\left(rac{Dp_{
m i}}{Dp_{
m i+1}}
ight)}$$

For fixed pore size tables (if selected), the following calculations are performed

(1) Average Fixed Pore Size (DF_{avg,j},A):

$$DF_{ ext{avg,j}} = rac{Dp_{F_{ ext{j}}} + D_{P_{F_{ ext{j}+1}}}}{2}$$

calculated for all intervals in the fixed pore size table.

For the intervals with between the Minimum BJH diameter and the Maximum BJH diameter.

(2) Cumulative Pore volume (VpF_{cum,i}, cm³/g):

$$V_P F_{ ext{cum.i}} = ext{INTERP}(D_P F_{ ext{i+1}})$$

where INTERP(x) is the value interpolated from the function $X=dP_{j+1}$ and $Y=VP_{cum,i}$, using an AKIMA semi-spline interpolation.

(3) Incremental Pore Volume (VpF_i, cum³/g):

$$V_P F_i = V_P F_{\mathrm{cum,i}} - V_P F_{\mathrm{cum,i-1}}$$

where $VpF_{cum,0}=0$



(4) Cumulative Surface Area (SAF_{cum,i}, m²/g):

$$SAF_{\mathrm{cum,i}} = \mathrm{INTERP}(D_P F_{\mathrm{i+1}})$$

where INTERP(x) is the value interpolated from the function $X = Dp_{j+1}$ and $Y = SA_{cum,j}$.

(5) Incremental Surface Area (SAF_i, m²/g):

$$SAF_{i} = SAF_{\text{cum,i}} - SAF_{\text{cum,i-1}}$$

where $SAF_{cum.0} = 0$.

(6) dV/dD pore volume ($dV/dDpF_i$, cm³/g-A):

$$rac{dV}{dDpF_{
m i}} = {
m INTERP}(DpF_{
m i+1})$$

where INTERP(x) is the value interpolated from the function $X = D_{avg,j}$ and $Y = dV/dD_i$

(7) dV/d log(D) pore volume [$dV/d log(DpF_i)$, cm³/g]:

$$rac{dV}{d\log(DpF_{
m i})} = {
m INTERP}(DpF_{
m i+1})$$

where INTERP(x) is the value interpolated from the function $X = D_{avg,j}$ and $Y = dV/d \log(D)_{i}$

(8) dA/dD pore area ($dA/dDpF_i$, m²/g-A):

$$rac{dA}{dDpF_{
m i}}={
m INTERP}(DpF_{
m i+1})$$

where INTERP(x) is the value interpolated from the function $X = D_{avg,j}$ and $Y = dA/dD_i$.

(9) dA/d log(D) pore area $[dA/d log(DpF_i), m^2/g]$:

$$rac{dA}{d\log(DpF_{
m i})} = {
m INTERP}(DpF_{
m i+1})$$

where INTERP(x) is the value interpolated from the function $X = D_{avg,j}$ and $Y = dA/d log(D)_i$.



COMPENDIUM OF VARIABLES

 ΔTd = thickness of layer desorbed from walls of newly opened pores (Å)

 ΔTw = thickness of adsorbed layer desorbed during interval (Å)

A = adsorbate property factor; from the BJH Adsorptive Options window

CSA = analysis gas molecular cross-sectional area (nm²), user-entered on the Adsorpt-

ive Properties window

CSAa = annular cross-sectional area of the desorbed layer (cm²)

CSAc = cross-sectional area of opening of newly opened pores (cm²)

 D_{avg} = average pore diameter (Å) Dp = pore (or core) diameter (Å)

F = fraction of pores open at both ends; from the BJH Adsorption Report Options win-

dow or the BJH Desorption Report Options window

LP = length of pore (cm/g)
P = relative pressure

Q = quantity adsorbed expressed as a volume (cm^3/g STP)

Rc = Kelvin radius (Å) of core

SAw = total surface area of walls exposed (cm²/g)
Tw = thickness of remaining adsorbed wall (Å)

Vc = volume desorbed from cores of newly opened pores (cm³/g)

Vd = volume of gas desorbed from walls of previously opened pores (cm³/g)

VI = liquid equivalent volume of volume adsorbed (cm³/g) V_{mol} = liquid molar volume, from the fluid property information



CRYSTALLITE SIZE FOR CHEMICAL ADSORPTION

$$d_{ ext{xtal}} = rac{1000k}{\overline{
ho}A_{ ext{metal}}}$$

Where

k = shape factor; 6 for sphere, 5 for cube

 $\overline{\rho}$ = weighted average density of the active metals



DATA SMOOTHING FOR CHEMICAL ADSORPTION

The data smoothing feature of the chemical adsorption program was developed to cope with noise and drift of pressure readings at very low pressures — usually at pressures well under 1 mmHg and very low gas volumes adsorbed. Data smoothing in the chemical adsorption program requires eleven points to satisfy the technique employed. However, most chemical adsorption analyses are performed with fewer than eleven points and at pressures higher than 1 mmHg, and so the data smoothing option is turned off by default. It may be enabled if required

TECHNIQUE

The analysis information consists of absolute pressure and volume adsorbed data pairs (P,V). In order to account for negative transducer shift, all coordinate pairs are scanned for points with pressures less than or equal to zero. If such points exist, then all pressure data are shifted by an amount large enough to convert the smallest pressure to a nominally small positive value.

After examination of the low pressure data is complete, the entire data set is subdivided into equally spaced intervals along the volume axis and corresponding pressures for these volumes are interpolated. The pressure data are smoothed by passing them through an eleven-point Blackman¹) convolution array. Once the smoothing is complete, the volume adsorbed values from the original data set are used to interpolate back to the corresponding smoothed analysis pressures. This amended set of data, consisting of the original volume adsorbed values and the fully smoothed pressure values, forms the isotherm data from which all subsequent report data are based. Coefficients for the Blackman convolution array are given below.

```
P_{\text{smoothed}} = 0.0029918P_{i-5} + 0.020005P_{i-4} + 0.0618580P_{i-3} + 0.1250300P_{i-2} + 0.1850900P_{i-1} + 0.2100500P_{i} + 0.1850900P_{i+1} + 0.1250300P_{i+2} + 0.061850P_{i+3} + (0.0200050)P_{i+4} + 0.0029918P_{i-5}
```

where

 P_{smoothed} = i^{th} smoothed pressure P_i pressure coordinate from the i^{th} equally spaced data point

¹⁾ Blackman, R.B. and Tukey, J.S., "The Measurement of Power Spectra," Dover (1958).



DFT (DENSITY FUNCTIONAL THEORY)

The adsorption isotherm is known to convey a great deal of information about the energetic heterogeneity and geometric topology of the sample under study. The data of physical adsorption have been used for many years as the basis for methods to characterize the surface area and porosity of adsorbents. Real solid surfaces rarely approach ideal uniformity of structure. It is accepted that in general, the surface of even a nonporous material presents areas of greater or lesser attraction for adsorbed molecules.

This energetic heterogeneity greatly affects the shape of the adsorption isotherm with the result that simple theories such as the Langmuir and BET formulas can, at best, give only approximate estimates of surface area. Porous solids virtually are never characterized by a single pore dimension, but instead exhibit a more or less wide distribution of sizes. The observed adsorption isotherm for a typical material is therefore the convolution of an adsorption process with the distribution of one or more properties which affect that process. This was first stated mathematically by Ross and Olivier 1) for the case of surface energy distribution and has become known as the integral equation of adsorption.

THE INTEGRAL EQUATION OF ADSORPTION

In a general form for a single component adsorptive, the integral equation of adsorption can be written as:

$$Q(p) = \int da \operatorname{dbdc...} q(p, a, b, c...) \int (a, b, c...)$$
(1)

where

Q(p) = the total quantity adsorbed per unit weight at pressure p,

a,b,c,... = a set of distributed properties,

f(a,b,c,...) = the distribution function of the properties, and

q(p,a,b,c,...) = the kernel function describing the adsorption isotherm on unit surface of

material with fixed properties a,b,c,...

Equation (1), a Fredholm integral of the first kind, is a member of a class of problems known as ill-posed, in that there are an infinite number of functional combinations inside the integral that will provide solutions. Even when the kernel function is known, experimental error in the data can make solving for even a single distribution function a difficult task. Solving for multiple distribution functions requires more data than provided by a single adsorption isotherm.

¹⁾ Ross and Olivier, J.P., "On Physical Adsorption," J. Wiley and Sons, New York (1964).



APPLICATION TO SURFACE ENERGY DISTRIBUTION

Under certain conditions, an energetically heterogeneous surface may be characterized by a distribution of adsorptive energies. The conditions are that the sample is not microporous, i.e., that adsorption is taking place on essentially a free surface with no pore filling processes at least to about 0.2 relative pressure. Secondly, that each energetically distinct patch contributes independently to the total adsorption isotherm in proportion to the fraction of the total surface that it represents. This condition is satisfied if the patches are relatively large compared to an adsorptive molecule, or if the energy gradient along the surface is not steep. In mathematical terms, this concept is expressed by the integral equation of adsorption in the following form:

$$Q(p) = \int d\epsilon \ q(p,\epsilon) \int (\epsilon)$$
 (2)

where

Q(p) = the experimental quantity adsorbed per gram at pressure p,

 $q(p,\varepsilon)$ = the quantity adsorbed per unit area at the same pressure, p, on an ideal free surface of energy ε , and

 $f(\varepsilon)$ = the total area of surface of energy ε in the sample.

The exact form of the energy-dependent term depends on the form of the model isotherms expressed in the kernel function and is provided in the model description.

APPLICATION TO PORE SIZE DISTRIBUTION

Similarly, a sample of porous material may be characterized by its distribution of pore sizes. It is assumed in this case that each pore acts independently. Each pore size present then contributes to the total adsorption isotherm in proportion to the fraction of the total area of the sample that it represents. Mathematically, this relation is expressed by:

$$Q(p) = \int dH \ q(p, H) \int (H) \tag{3}$$

where

Q(p) = the experimental quantity adsorbed at pressure p,

q(p,H) = the quantity adsorbed per unit area at the same pressure, p, in an ideal pore of size H, and

f(H) = the total area of pores of size H in the sample.



Numerical values for the kernel functions in the form of model isotherms can be derived from modern statistical mechanics such as density functional theory or molecular simulations, or can be calculated from one of various classical theories based on the Kelvin equation. Several types are found in the models library.

PERFORMING THE DECONVOLUTION

The integrations in equations (2) and (3) are carried out over all surface energies or pore sizes in the model. The functions $q(p,\varepsilon)$ and q(p,H), which we call the kernel functions, are contained in numeric form as model isotherms. Because, in general, there is no analytic solution for equation (1), the problem is best solved in a discrete form; the integral equation for any distributed property Z becomes a summation:

$$Q(p) = \sum_{i} q(p, Z_i) \int (Z_i)$$
(4)

Given a set of model isotherms, q(p,Z), from a model chosen from the models library and an experimental isotherm, Q(p), contained in a sample information file, the software determines the set of positive values f(Z) that most nearly, in a least squares sense, solves equation (4). The distributed property, surface energy or pore size, is then displayed on the *Report Options* window as a selection of tables or graphs.

REGULARIZATION

DFT allows a selectable regularization (also referred to as smoothing) constraint to be applied during the deconvolution process to avoid over-fitting in the case of noisy data or ill-fitting models. The method used is based on co-minimization of the second derivative of the distribution. The relative weight given to this term is determined by the value of the regularization parameter, which is set on the *DFT Pore Size* or *Surface Energy* window and also is shown in the header of reports. The value of the regularization parameter varies from zero (for no second derivative constraint) to ten (indicating a weight equal to minimizing the residuals), or even larger. When the distribution and residuals obtained change little with the value of the regularization parameter, it indicates that the chosen model provides a good representation of the data. Conversely, a large sensitivity to the regularization parameter might indicate inadequate data or a poor choice of model to represent the data.



DOLLIMORE-HEAL PORE VOLUME AND AREA DISTRIBUTION

The calculations for the Dollimore-Heal reports are the same as those for BJH, except for the calculation of average pore diameter and pore length.

PORE DIAMETER

Pore diameter is determined from the Kelvin radius and thickness equation:

$$D_{\mathrm{i}} = 2r_{\mathrm{k}}P_{i} + tP_{\mathrm{i}}$$

The average pore diameter is the arithmetic mean of the diameters that bound the interval.

$$\overline{D_{
m i}} = \left(rac{D_i + D_{
m i+1}}{2}
ight)$$

PORE LENGTH

$$l_{
m i}=rac{A_{
m p,i}+10^8}{\pi\overline{D}_{
m i}}$$

$$A_{
m p,i} = rac{4 imes (10^8 \Delta V_{
m p})}{\overline{D_{
m i}}}$$

$$\Delta V_{
m p} = C
u (D(Q_{
m i-1}-Qi) - \Delta t imes 10^8 (A_{
m p,cum} - 2\pi ar{t} l_{
m i,cum}))$$

$$Cv=\left(rac{ar{D_{\mathrm{i}}}}{2(ar{\pi}_{\mathrm{k}}+t(P_{\mathrm{i}})-t(P_{\mathrm{i+1}}))}
ight)^{2}$$

$$ar{t}=rac{ar{D_{
m i}}}{2-ar{r}_{
m k}}$$

$$ar{r}_{ ext{k}} = rac{(r_{ ext{k,i}} + r_{ ext{k,i+1}})}{2}$$



where

 ΔV_p = Change in pore volume

 $A_{p,i}$ = Pore surface area

 $A_{p,i,cum}$, $I_{i,cum}$ = Summations over the lengths and areas calculated so far

Cv = Volume correction factor D = Density conversion factor $ar{r}_k$ = Average Kelvin radius $ar{t}$ = Average thickness



DUBININ-ASTAKHOV

The Dubinin-Astakhov equation is:

$$\log(Q) = \log(Q_0) - \left(rac{RT}{eta E_0}
ight)^N imes \left(\lograc{P_0}{P}
ight)^N$$

where

β = the affinity coefficient of the analysis gas relative to the P₀ gas, from the *Dubinin*Adsorptive Options window

 E_0 = characteristic energy (*kj/mol*)

N = Astakhov exponent, may be optimized or user entered from the *Dubinin Report Options* window

P = equilibrium pressure

 P_0 = saturation vapor pressure of gas at temperature T

Q = quantity adsorbed at equilibrium pressure (cm³/g STP)

 Q_0 = the micropore capacity (cm³/g STP)

R = the gas constant (0.0083144 kj/mol)

T = analysis bath temperature (K)

For each point designated for Dubinin-Astakhov calculations, the following calculations are done:

$$LV = \log(Q)$$

$$LP = \log(rac{P_0}{P})^N$$

A least-squares fit is performed on the (LP,LV) designated pairs where LP is the independent variable and LV is the dependent variable. If the user selected Yes for the Optimize Astakhov Exponent prompt, a systematic search for the optimum value of N is conducted by recalculating the linear regression and selecting the value of N that gives the smallest standard error of the y-intercept. The exponent N is optimized to within 10^{-4} . If the optimum value for N is not found in this range, an exponent of 2 is used. The following are calculated:



- a. Slope (S cm³/g STP)
- b. Y-intercept (YI cm³/g STP)
- c. Error of the slope ($S_{err} cm^3/g STP$)
- d. Error of the y-intercept (YI_{err} cm 3 /g STP)
- e. Correlation coefficient
- f. Optimized Astakhov exponent (N)

Using the results of the above calculations, the following can be calculated:

Monolayer Capacity (cm³/g STP):

$$Q_0 = 10^{YI}$$

Micropore Volume (cm³/g):

$$V_{
m i}=rac{Q_i V_{
m mol}}{22414}$$

where

 V_{mol} = liquid molar volume conversion factor from the fluid property information

Limiting Micropore Volume (cm³/g):

$$V_0 = \frac{Q_0 V_{
m mol}}{22414 {
m cm}^3 {
m STP}}$$

where

 V_{mol} = liquid molar volume from the fluid property information

Error of Limiting Micropore Volume (cm³/g):

$$V_{0,\mathrm{err}} = W_0 (10 Y I_{\mathrm{err}} - 1.0)$$

Characteristic Energy (KJ/mol):

$$E = rac{2.303(RT)}{eta(2.303 imes S)^{1/{
m N}}}$$



Modal Equivalent Pore Diameter (Å):

$$D_{\mathrm{mode}} = 2 \Bigg\{ \left[rac{3N}{3N+1}
ight]^{1/N} imes \left[rac{10^3 n m^3 / {\dot A}^3}{eta \cdot E_0}
ight] \Bigg\}^{1/3}$$

where

β = affinity coefficient of the analysis gas relative to the P₀ gas from the *Dubinin Adsorptive Options* window

Maximum Differential Pore Volume (cm³/g-Å):

This value is also known as frequency of the mode¹).

$$rac{dV}{dD_{
m mode}}{
m Max} = 0.5(3N+1)W_0igg[rac{3N+1}{3N}igg]^{1/3N}igg[rac{eta\cdot E_0}{((10^3nm^3)/{\dot A}^3)}igg]^{1/3} {
m exp}igg(-igg[rac{3N+1}{3N}igg]igg)$$

Mean Equivalent Pore Width (Å):

$$D_{
m mean} = 2 imes rac{\left[rac{(10^3 {
m nm}^3) \dot{A}^3}{eta \cdot E_0}
ight]^{1/3}}{ au(rac{3N+1}{3N})}$$

Micropore surface area (m²/g):

$$SDA = 1000 imes 2.0 imes W_0 imes \left[rac{E_0}{k}
ight]^{1/3} imes au \left(rac{3N+1}{3N}
ight)$$

 Γ is calculated by a polynomial approximation over the domain $0 \le x \le 1$:

$$au(x+1) = 1 + b_1 x + b_2 x^2 + b_3 x^3 + b_4 x^4 + b_5 x^5 + b_6 x^6 + b_7 x^7 + b_8 x^8 \ + arepsilon \mathbf{x} |arepsilon \mathbf{x}| \leq 3 (10^{-7})$$

¹⁾ Ross and Olivier, J.P., "On Physical Adsorption," J. Wiley and Sons, New York (1964).



where

 b_1 -0.57719 1652 b_2 0.98820 5891 b_3 -0.89705 6937 0.91820 6857 b₄ -0.75670 4078 b_5 0.48219 9394 b_6 -0.19352 7818 b_7 0.03586 8343 bg

and where

$$x+1=\left(rac{3N+1}{3N}
ight)$$

Equivalent Pore Diameter (Å):

$$D_{
m i} = 2 igg[rac{-ig(rac{10^3 {
m mm}^3/{
m \AA}^3}{eta \cdot E_0}ig)^N}{{
m ln}(W_{
m i}) - {
m ln}(W_0)} igg]^{1/3N}$$

dV/dD Pore Volume (cm³/g-Å):

$$egin{aligned} rac{dV}{dD_{\mathrm{i}}} &= 0.5 imes W_0 imes 3Nigg(rac{10^3 \mathrm{nm}^3/\mathrm{\dot{A}}^{3N}}{eta \cdot E_0}igg)igg(rac{D_{\mathrm{i}}}{2}igg)^{-(3N+1)} \ & imes \expigg[-igg(rac{10^3 \mathrm{nm}^3/\mathrm{\dot{A}}^3}{eta \cdot E_0}igg)^Nigg(rac{D_{\mathrm{i}}}{2}igg)^{-3N}igg] \end{aligned}$$



DUBININ-RADUSHKEVICH

The Dubinin-Radushkevich¹) equation is:

$$\log(Q) = \log(Q_0) - rac{BT^2}{eta} imes \left[\lograc{P_0}{P}
ight]^2$$

where

 β = the affinity coefficient of analysis gas relative to P₀ gas (for this application β is taken to be 1)

to be 1)

B = a constant

 P_0 = saturation vapor pressure of gas at temperature T

P = equilibrium pressure

Q = quantity adsorbed at equilibrium pressure (cm³/g STP)

 Q_0 = the micropore capacity (cm³/g STP)

T = analysis bath temperature (K), from the P_0 and T emperature Options window

For each point designated for Dubinin-Radushkevich calculations, the following calculations are done:

$$LV = \log(Q)$$

$$LP = \log(rac{P_0}{P})^2$$

The intercept, log(Vo) can be found by performing a least-squares fit on the (LP,LV) designated pairs where LP is the independent variable and LV is the dependent variable. Assuming the adsorption of gas is restricted to a monolayer, Vo is the monolayer capacity. Based on this assumption, the following are calculated:

¹⁾ Dubinin, M., *Carbon 21*, 359 (1983); Dubinin, M., *Progress in Surface and Membrane Science 9*, 1, Academic Press, New York (1975); Dubinin, M. and Astakhov, V., *Adv. Chem. Ser. 102*, 69 (1971); Lamond, T. and Marsh, H., *Carbon 1*, 281, 293 (1964); Medek, J., *Fuel 56*, 131 (1977); Polanyi, M., *Trans. Faraday Soc. 28*, 316 (1932); Radushkevich, L., *Zh. fiz. Kemi. 33*, 2202 (1949); Stoeckli, H., *et al*, *Carbon 27*, 125 (1989).



- a. Slope (S cm³/g STP)
- b. Y-intercept (Y/cm^{3/}g STP)
- c. Error of the slope (S_{err} cm³/g STP)
- d. Error of the y-intercept (YI_{err} cm³/g STP)
- e. Correlation coefficient

Using the results of the above calculations, the following can be calculated:

Monolayer Capacity (cm³/g STP):

$$Q_0=10^{YI}$$

Error of Monolayer Capacity (cm³/g STP):

$$Q_{0,
m err} = Q_0 (10^{YI,
m err} - 1.0)$$

Micropore surface area (m²/g):

$$SDP = rac{\sigma Q_0 N_A}{22414 {
m cm}^3 \left(rac{10^{18} {
m mm}^2}{{
m m}^2}
ight)}$$

where

 σ = molecular cross sectional area of gas (nm²) from the *Adsorptive Properties* window



EQUATION OF STATE

The ideal gas law relates pressure, volume, temperature, and quantity of gas.

$$n=rac{PV}{RT}$$

where

P = pressure

V = volume

R = a constant that depends on the units of n

For
$$n \text{ in } cm^3$$
, STP, R = P_{STD} / T_{STD}

For *n* in moles, $R = 8.3145 \text{ J mol}^{-1} \text{ K}^{-1}$

T = temperature

The non-ideality correction is made by replacing P with $P(1 + \alpha P)$, where α is the non-ideality factor. This gives

$$n=rac{PV}{RT}(1+lpha P)$$

The real gas equation of state uses

$$n=rac{PV}{RT_z(P,T)}$$

where z(P,T) is the compressibility factor for the gas at the given pressure and temperature.



EQUILIBRATION

Equilibration is reached when the pressure change per equilibration time interval (first derivative) is less than 0.01% of the average pressure during the interval. Both the first derivative and average pressure are calculated using the Savitzky-Golay¹) convolution method for polynomial functions. The following equations are those used to compute weighted average and first derivative, respectively, for the 6th point of an 11-point window:

$$P_{\text{avg}} = \frac{-36(P_{11} + P_1) + 9(P_{10} + P_2) + 44(P_9 + P_3) + 69(P_8 + P_4) + 84(P_7 + P_5) + 89(P_6)}{429}$$

$$P_{
m chg} = rac{5(P_{11}-P_1)+4(P_{10}-P_2)+3(P_9-P_3)+2(P_8-P_4)+(P_7-P_5)}{110}$$

$$P_{\mathrm{pcp},i} = 100\% rac{P_{\mathrm{chg}}}{P_{\mathrm{avg}}}$$

pressure change per equilibration time interval

where the numerical constants are from the Savitzky-Golay convolution arrays, and

 P_{avg} = average pressure P_{chq} = change in pressure

P_{pcp,i} = percent change per interval

 P_i = ith pressure reading taken at equilibrium intervals



If a non-zero value that is too small is entered for the maximum equilibration time, the points are collected before equilibration is reached.



If P_{avg} is greater than 0.995 times the current P_0 , equilibration will not take place until the *Minimum equilibration delay for P/P₀ 0.995* has expired, in addition to the standard equilibration criteria.

¹⁾ Savitzky, A. and Golay, M.J.E., Anal. Chem. 36, 1627 (1964).



F-RATIO METHOD

The *f*-Ratio is the quantity adsorbed divided by the quantity adsorbed in a reference isotherm at the same pressure.

$$f_i = rac{Q_i}{Q_{ ext{ref}}P_i}$$

The reference quantity adsorbed is found by spline interpolation of the reference isotherm.



FREE SPACE FOR CHEMICAL ADSORPTION

$$V_{ ext{FS}} = rac{V_{ ext{LOW}}}{T_1} imes \left[rac{P_1}{P_2} - 1
ight] imes T_{ ext{STD}}$$

where

 P_1 = system manifold pressure before dosing helium onto sample (μ mHg)

 P_2 = system manifold pressure after dosing helium onto sample (μ mHg)

 T_1 = system manifold temperature before dosing helium onto sample (K)

 T_2 = system manifold temperature after raising Dewar and equilibrating with helium (K)

 T_{STD} = standard temperature (273.13 K)

 V_{FS} = volume of free space (cm³ at standard temperature)

 $V_{\rm LOW}$ = lower manifold volume (cm³)



FREUNDLICH ISOTHERM

The Freundlich isotherm has the form:

$$rac{Q}{Q_{\mathrm{m}}}=CP^{rac{1}{\mathrm{m}}}$$

where

C = temperature-dependent constant

m = temperature-dependent constant

P = equilibrated collected pressure measured by gauge at temp T_{amb}

Q = quantity of gas adsorbed

 Q_m = quantity of gas in a monolayer

The pressure is absolute; typically, m > 1. In terms of quantity adsorbed,

$$Q=Q_{
m m}CP^{rac{1}{
m m}}$$

Taking the log of both sides yields

$$\log Q = \log Q_{\mathrm{m}}C + \frac{1}{\mathrm{m}}\log P$$



HEAT OF ADSORPTION FOR PHYSICAL ADSORPTION

The adsorption isostere is represented by

$$\ln\left(rac{P}{P_0}
ight) = rac{q_i}{RT} + C$$

where

 q_i = isosteric heat of adsorption

C = unknown constant

R = gas constant

T = analysis temperature

The isosteric heat of adsorption, q_i , is determined by finding the slope of $ln (P/P_0)$ as a function of 1/RT for a set of isotherms measured at different temperatures.



HORVATH-KAWAZOE

A relative pressure lower limit is determined such that L- d_0 never equals zero. All pressure points less than this limit are discarded. For each collected relative pressure point, values of L are chosen in an iterative manner, and the relative pressure (P/P_0) determined by solving one of the following equations:

- Slit Pore Geometry (original Horvath-Kawazoe)
- Cylinder Pore Geometry (Saito / Foley)
- Sphere Pore Geometry (Cheng / Yang)

SLIT PORE GEOMETRY (ORIGINAL HORVATH-KAWAZOE)

When using the original Horvath-Kawazoe¹) method, the following equation is solved for each value of P. The value of L is determined when the solved-for relative pressure is within 0.1% of the collected absolute pressure:

$$\ln rac{P}{P_0} = rac{N_A}{RT} rac{IP imes 10^{32}}{\sigma^4 L - 2d_0} rac{\sigma_4}{3L - {d_0}^3} - rac{\sigma^{10}}{9L - {d_0}^9} - rac{\sigma^4}{3{d_0}^3} + rac{\sigma^{10}}{9{d_0}^9}$$

where

 10^{32} = the number of cm⁴ that are equal to Å⁴

 σ = gas solid nuclear separation at zero interaction energy (Å), $\frac{Z_s + Z_A}{2}$

 $d_0 = \frac{D_A + D_s}{2}$

where

D_A = molecular diameter (Å) from the *Horvath-Kawazoe Physical Properties* window

D_S = diameter of sample atom (Å) from the *Horvath-Kawazoe Physical Properties* window

IP = interaction parameter (erg-cm⁴) from the *Horvath-Kawazoe Report Options* window

 N_A = Avogadro Constant (6.02214129 × 10²³ mol⁻¹)

L = pore width (nucleus to nucleus) (Å)

P = equilibrium pressure

¹⁾ Horvath, G. and Kawazoe, K., J. Chem. Eng. Japan 16(6), 470 (1983).



 P_0 = saturation pressure

R = gas constant $(8.31441 \times 10^7 \text{ erg/mol } K)$

T = analysis bath temperature (K), from an entered or calculated value on the P_0 and Temperature Options window

where:

Z_s = sample equilibrium diameter at zero interaction energy (Å) from the *Horvath-Kawazoe Physical Properties* window

Z_A = zero interaction energy diameter from the *Horvath-Kawazoe Physical Properties* window

CYLINDER PORE GEOMETRY (SAITO/FOLEY)

When using the Saito/Foley¹) method, the following equation is solved for each value of P. The value of L is determined when the solved-for relative pressure is within 0.1% of the collected absolute pressure:

$$egin{split} \ln\left(rac{P}{P_0}
ight) &= rac{3}{4}rac{\pi N_A}{RT} imes rac{IP imes 10^{32}}{{d_0}^4} imes \sum_{k=0}^{\infty} \left[rac{1}{\mathrm{k}+1}igg(1-rac{d_0}{r_\mathrm{p}}igg)^{2\mathrm{k}} imes
ight. \ \left. \left\{rac{21}{32}lpha_\mathrm{k}igg(rac{d_0}{r_\mathrm{p}}igg)^{10} - eta_\mathrm{k}igg(rac{d_0}{r_\mathrm{p}}igg)^4
ight\}
ight] \end{split}$$

where

 10^{32} = the number of Å⁴ that are equal to 1 cm⁴

$$\beta_k = \left(\frac{-1.5-k}{k}\right)^2 \beta_{k-1}, \beta_0 = 1.0$$

$$\alpha_{k}$$
 = $\left(\frac{-4.5-k}{k}\right)^{2}\alpha_{k-1}, \alpha_{0}=1.0$

¹⁾ Saito, A. and Foley, H. C., AIChE Journal 37(3), 429 (1991).



$$d_0 = \frac{D_A D_s}{2}$$

where

D_A = molecular diameter (Å) from the *Horvath-Kawazoe Physical Properties* window

D_S = diameter of sample atom (Å) from the *Horvath-Kawazoe Physical Properties* window

IP = interaction parameter (10⁻⁴³ erg-cm⁴) from the *Horvath-Kawazoe Report Options* window

 N_A = Avogadro Constant (6.02214129 × 10²³ mol⁻¹)

L = pore width (nucleus to nucleus) (Å)

P = equilibrium pressure

P_o = saturation pressure

R = gas constant $(8.31441 \times 10^7 \text{ erg/mol } K)$

 r_p = radius of the cylindrical pore, $\frac{L}{2}$

T = analysis bath temperature (K), from an entered or calculated value on the Po and Temperature Options window



SPHERE PORE GEOMETRY (CHENG/YANG)

When using the Cheng / Yang 1) method, the following equation is solved for each value of P. The value of L is determined when the solved-for relative pressure is within 0.1% of the collected absolute pressure:

$$\ln\left(rac{P}{P_0}
ight) = rac{6N_1\epsilon^*_{12} + N_2\epsilon^*_{22}L^3 imes 10^{32}}{RTL - d_0{}^3} igg[-\left(rac{d_0}{L}
ight)^6 igg(rac{T_1}{12} + rac{T_2}{8}igg) + \left(rac{d_0}{L}
ight)^{12} igg(rac{T_3}{90} + rac{T_4}{80}igg) igg]$$

where

 10^{32} = the number of cm⁴ that are equal to Å⁴

$$\epsilon^*_{12} = \frac{\dot{A}_S}{4d_s^6}, \text{where} \dot{A}_s = \frac{6(mc^2)\alpha_S\alpha_A}{\frac{\alpha_S}{X_S} + \frac{\alpha_A}{X_A}}$$

$$\epsilon^*_{22} = \frac{\dot{A}_A}{4D_A^6}$$
, where $\dot{A}_A = \frac{3(mc^2)(\alpha_A)(\chi_A)}{2}$

$$d_0 = \frac{D_A + D_S}{2}$$

where

D_A = molecular diameter (Å) from the *Horvath-Kawazoe Physical Properties* window

D_S = diameter of sample atom (Å) from the *Horvath-Kawazoe Physical Properties* window

L = pore width (nucleus to nucleus) (Å)

 $N_1 = 4\pi L_2 N_S$, where $N_s =$ number of sample atoms/cm² at monolayer

 $N_2 = 4\pi (L-d_0)^2 N_A$, where $N_S = \text{number of gas molecules/cm}^2$

P = equilibrium pressure

 P_0 = saturation pressure

R = gas constant $(8.31441 \times 10^7 \text{ erg/mol K})$

T = $\frac{\text{analysis bath temperature } (K), \text{ from an entered or calculated value on the } P_0 \text{ and } Temperature Options}{\text{Notion of the } P_0 \text{ and } P_0 \text$

¹⁾ Cheng, Linda S. and Yang, Ralph T., Chemical Engineering Science 49(16), 2599-2609 (1994).



$$T_1 = rac{1}{(1-S)^3} - rac{1}{(1+S)^3}$$

$$T_2 = rac{1}{(1+S)^2} - rac{1}{(1-S)^2}$$

$$T_3 = rac{1}{(1-S)^9} - rac{1}{(1+S)^9}$$

$$T_4 = rac{1}{(1+S)^8} - rac{1}{(1-S)^8}$$

where
$$S=rac{L-d_0}{L}$$

CHENG/YANG CORRECTION

This factor corrects for the nonlinearity of the isotherm. It adds an additional term to the equations for the different geometrics:

$$\ln\left(rac{P}{P_0}
ight) = G(L) - \left[1 - rac{1}{ heta} \mathrm{ln}\left(rac{1}{1 - heta}
ight)
ight]$$

where

G(L) = one of the Horvath-Kawazoe equations given above

 θ = degree of void filling; θ is estimated by first computing the monolayer capacity (Q_m) with the Langmuir equation over the range of data points from relative pressure 0.02 to 0.2 or the maximum relative pressure included in the Horvath-Kawazoe analysis. θ is computed as the quantity adsorbed over Q_m .



INTERACTION PARAMETER

The interaction parameter (IP) results from the following calculations:

The Kirkwood-Muller dispersion coefficients

$$A_{
m S}=rac{6mc^2lpha_{
m S}lpha_{
m A}}{rac{lpha_{
m S}}{\chi_{
m S}}+rac{lpha_{
m A}}{\chi_{
m A}}}$$

$$A_{
m A}=rac{3mc^2lpha_{
m A}\chi_{
m A}}{2}$$

where

 α_A = polarizability of gas molecule (cm³)

 $\alpha_{\rm S}$ = polarizability of sample atoms (cm³)

 mc^2 = kinetic energy of electron (0.8183 × 10⁻⁶ erg)

 X_A = diamagnetic susceptibility of gas molecule (cm³)

$$IP = (N_{
m M}A_{
m A}) + (N_{
m S}A_{
m S})$$

where:

N_M = number of gas molecules/cm² at monolayer from the *Horvath-Kawazoe* Physical Properties window

N_S = number of sample atoms/cm² from the *Horvath-Kawazoe Physical Properties* window

 X_s = diamagnetic susceptibility of sample atom (cm³)

See Interaction Parameter Components Table on page 38 for recommended values.



ADDITIONAL CALCULATIONS

Based on the previous calculations, the following can be calculated:

Adjusted Pore Width (Å):

(Shell to Shell)

$$AL_{\rm i} = L_{\rm i} - D_{\rm S}$$

Cumulative Pore Volume (cm³/g):

$$V_{ ext{cum,i}} = rac{Q_{ ext{i}}V_{ ext{mol}}}{22414 ext{cm}^3 ext{STP}}$$

where

 V_{mol} = liquid molar volume from the fluid property information

dV/dD Pore Volume (cm³/g-Å):

$$rac{dV}{dD_{
m i}} = rac{V_{
m cum,i} - V_{
m cum,i-1}}{AL_{
m i} - AL_{
m i-1}}$$

Median Pore Width (A):

$$V_{
m half} = rac{V_{
m cum,n}}{2}$$

$$D_{
m med} = \expigg(\ln D_{
m l} + (\ln D_{
m g} - \ln D_{
m l}) rac{\ln V_{
m half} - \ln V_{
m l}}{\ln V_{
m g} - \ln V_{
m l}}igg)$$

where

 D_{l} = pore width (L_{i}) that corresponds to V_{l}

 D_g = pore width (L_i) that corresponds to V_g

 $V_{\text{cum,n}}$ = total cumulative pore volume ($V_{\text{cum,i}}$) for points designated for Horvath-Kawazoe

calculations

 V_q = cumulative pore volume ($V_{\text{cum,i}}$) for first point greater than V_{half}

 V_{half} = 50% of total cumulative pore volume

 V_I = cumulative pore volume ($V_{\text{cum,i}}$) for first point less than V_{half}



Interaction Parameter Components Table

Gas	Bath Temperature (K)	Sample Type	Interaction Parameter Calculated Value *
Argon	87.3	Carbon (Ross/Olivier value) Carbon (Horvath/Kawazoe value) Zeolite	2.61 5.89 3.19
Carbon Dioxide	298.15	Carbon (Ross/Olivier value) Carbon (Horvath/Kawazoe value) Zeolite	4.20 9.20 5.08
	273.15	Carbon (Ross/Olivier value) Carbon (Horvath/Kawazoe value) Zeolite	4.34 9.35 5.22
	194.65	Carbon (Ross/Olivier value) Carbon (Horvath/Kawazoe value) Zeolite	4.72 9.72 5.60
Nitrogen	77.15	Carbon (Ross/Olivier value) Carbon (Horvath/Kawazoe value) Zeolite	2.84 6.53 3.49

^{*} The interaction parameter is entered in the *Horvath-Kawazoe Report Options* window in the following field:

Interaction parameter: (calculated value) \times 10⁻⁴³ erg-cm⁴



The following values were used to calculate the values in the previous table.

Carbon-Graphite

D_S = 3.40 N_S = 3.845 × 10¹⁵ X_S = 1.05 × 10⁻²⁹ (Ross/Olivier) 13.5 × 10⁻²⁹ (Horvath/Kawazoe, implicit) α_S = 1.02 × 10⁻²⁴

Zeolite

$$D_S$$
 = 3.04
 N_S = 3.75 × 10¹⁵
 X_S = 1.94 × 10⁻²⁹
 α_S = 0.85 × 10⁻²⁴

Nitrogen

$$D_A$$
 = 3.00
 N_A = 6.71 × 10¹⁴
 X_A = 3.6 × 10⁻²⁹
 α_A = 1.76 × 10⁻²⁴

Argon

$$D_A$$
 = 2.95
 N_A = 7.608 × 10¹⁴
 X_A = 3.22 × 10⁻²⁹
 α_A = 1.63 × 10⁻²⁴

Carbon Dioxide

$$D_A = 3.23$$

$$N_A = 4.567 \times 10^{14} (25 \,^{\circ}\text{C})$$

$$5.45 \times 10^{14} (0 \,^{\circ}\text{C})$$

$$7.697 \times 10^{14} (-78 \,^{\circ}\text{C})$$

$$X_A = 5.0 \times 10^{-29}$$

$$\alpha_A = 2.7 \times 10^{-24}$$

 D_A values are from van der Waal's constant.

 N_A values are from liquid densities.

X and α values are derived from data found in Ross and Olivier¹).

¹⁾ Ross and Olivier, J.P., "On Physical Adsorption," J. Wiley and Sons, New York (1964)



The physical parameters referenced in Saito/Foley are:

Aluminophosphate

$$D_S$$
 = 2.60
 N_S = 1.48 × 10¹⁵
 X_S = 1.3 × 10⁻²⁹
 α_s = 2.5 × 10⁻²⁴

Aluminosilicate

$$D_S$$
 = 2.76
 N_S = 1.31 × 10¹⁵
 X_S = 1.3 × 10⁻²⁹
 α_S = 2.5 × 10⁻²⁴



LANGMUIR SURFACE AREA FOR CHEMICAL ADSORPTION

TRANSFORM FOR CHEMICAL ADSORPTION

The Langmuir isotherm is

$$\frac{Q}{Q_m} = \frac{bP}{1+bP}$$

The isotherm is transformed so that P/Q is plotted as a function of pressure. The transformed data are fitted with a straight line. the slope (m) and intercept (y_0) of the fit line are used in the calculations below.

SURFACE AREA

$$A_{ ext{Lang}} = rac{ar{A}_{ ext{atom}}ar{S}N_A}{V_{ ext{mol}}m} \cdot 10^{-18} rac{m^2}{nm^2}$$

MONOLAYER CAPACITY

$$Q_m = \frac{1}{m}$$

LANGMUIR b VALUE

$$b=rac{1}{y_0Q_m}$$

DISSOCIATIVE CHEMICAL ADSORPTION

The Langmuir isotherm may be derived for dissociative chemical adsorption.

$$rac{Q}{Q_m}=rac{b\sqrt{P}}{1+b\sqrt{P}}$$

The calculations are performed with the slope and intercept of a fit of $\frac{b\sqrt{P}}{Q}$ as a function of \sqrt{P} .



LANGMUIR SURFACE AREA FOR PHYSICAL ADSORPTION

For each point designated for surface area calculations, the Langmuir 1) transformation is calculated as:

$$L=rac{P_{
m rel}}{N_{
m ads}}$$

where L is in units of g/cm³ STP.

A least-squares fit is performed on the (P_{rel}, L) designated pairs where P_{rel} is the independent variable and L is the dependent variable. The following are calculated:

- a. Slope (S g/cm³ STP)
- b. Y-intercept (Y_{int} g/cm³ STP)
- c. Error of the slope (S_{err} g/cm³ STP)
- d. Error of the y-intercept (YI_{err} g/cm³ STP)
- e. Correlation coefficient

Using the results of the above calculations, the following can be calculated:

Langmuir Surface Area (m²/g):

$$SA_{
m LAN} = rac{{
m CSA_{ imes}N_A}}{(22414{
m cm}^3{
m STP})(rac{10^{18}nm^2}{m^2})S}$$

where

CSA = analysis gas molecular cross-sectional area (nm²), user-entered on the Adsorptive Properties window

 N_A = Avogadro constant (6.02214129 × 10²³ mol⁻¹)

Quantity of the Monolayer (cm³/g STP):

$$Q_m = \frac{1}{s}$$

¹⁾ Langmuir, I., J. *Am. Chem. Soc. 38*, 2267 (1916); J. Am. Chem. Soc. 40, 1361 (1918); Phys. Rev 8, 149 (1916).



Langmuir b Value:

$$b = Y_{\mathrm{int}} V_m$$

Error of the Langmuir Surface Area (m²/g):

$$LAN_{
m err}=rac{SA_{
m Lan}S_{
m err}}{S}$$



METAL DISPERSION FOR CHEMICAL ADSORPTION

The metal dispersion is the percentage of the active metal available for interaction with the adsorbate.

$$\%M_{
m DISP} = rac{100\% imes100\%}{22414^*} imes rac{Q_0 imes SF_{
m CALC}}{rac{W_{
m atomic1}}{W_{
m ATOMIC2}} + rac{W_{
m weight3}}{W_{
m ATOMIC3}} + rac{W_{
m weight3}}{W_{
m ATOMIC3}}}$$

where

 $%M_{\text{DISP}}$ = metal dispersion (%)

% weight = % of sample weight for first metal

% weight = % of sample weight for second metal

%weight₃ = % of sample weight for third metal

 Q_0 = quantity adsorbed intercept of the best fit to the primary, repeat, or dif-

ference data. For the Sinfelt report, Q₀ is the difference of the primary and

repeat intercepts.

 SF_{CALC} = calculated stoichiometry factor

 W_{ATOMIC1} = atomic weight of first metal (g/mole)

 W_{ATOMIC2} = atomic weight of second metal (g/mole)

 W_{ATOMIC3} = atomic weight of third metal (g/mole)

^{*} The volume one mole of gas occupies (cm³ STP/mole of gas).



METALLIC SURFACE AREA

The metallic surface area per gram of sample is the total active metal surface area available for interaction with the adsorbate.

$$ext{Msa} = rac{6.023 imes10^{23}}{22414^*} imes V imes SF_{ ext{CALC}} imes A_{ ext{AREA}}$$

where

 $A_{\Delta R F \Delta}$ = effective area of 1 active metal atom (m²/atom)

Msa = metallic surface area (m²/g sample)

 SF_{CALC} = calculated stoichiometry factor

 $V = (cm^3/g STP)$

If Msa is being calculated from analysis data, then V is the volume intercept derived form the best line fit of the points selected for line fit by the user. If Msa is being calculated from difference data, then V is the average of the volume

differences (VDIFF) for the points selected for line fit.

 V_{DIFF} = difference in volume between the first analysis and the repeat analysis

Assuming that all species of metals disperse evenly, the metallic surface area per gram of metal is

$$N_{
m sa} = MSAigg(rac{100\%}{\% {
m weight1} + \% {
m weight2} + \% {
m weight3}}igg)$$

where

%weight1 = % of sample weight for first metal
 %weight2 = % of sample weight for second metal
 %weight3 = % of sample weight for third metal
 MSA = metallic surface area (m²/g sample)
 Nsa = metallic surface area (m²/g metal)

^{*} The volume one mole of gas occupies (cm³ STP/mole of gas)



MP-METHOD

With the (t_i,Q_i) data pairs 1), the Akima semi-spline interpolation method is used to interpolate quantity adsorbed values based on thickness values that are evenly spaced 0.2 angstrom apart starting at the first outlier point. Outliers are defined as those points that have the maximum instantaneous slope within an iteratively shrinking subset of all points. The remaining pore surface area calculation result is the slope of the line defined by two consecutive interpolated points. The slopes of each pair of consecutive points from the origin to the last point must be monotonically decreasing and non-negative. With the interpolated points set the following can be calculated:

Average pore hydraulic radius (Å):

$$R_{
m i}=rac{t_{
m i}+t_{
m i-1}}{2}$$

Remaining pore surface area for the ith point (m²/g):

$$S_{
m i} = rac{Q_{
m i} - Q_{
m i-1}}{t_{
m i} - t_{
m i-1}} rac{V_{
m mol}}{22414 {
m cm}^3 {
m STP}} imes 10^4$$

where

 10^4 = unit conversions

 V_{mol} = liquid molar volume from the fluid property information

Incremental pore surface area occluded for the i^{th} point (m²/g):

$$S_{\mathrm{inc},i} = S_{\mathrm{i-1}} - S_{\mathrm{i}}$$

Cumulative pore surface area occluded for the ith point (m²/g):

$$S_{\text{cum}_i} = S_{\text{inc,i}} + S_{\text{inc,i-1}} + \ldots + S_{\text{inc,i}}$$

dA/dR pore surface area for the ith point (m²/g-Å):

$$rac{dA}{dR_{
m i}}=rac{S_{
m inc_i}}{t_{
m i}-t_{
m i-1}}$$

¹⁾ Mikhail, R., Brunauer, S. and Bodor, E., J. Colloid and Interface Sci. 24, 45 (1968).



Incremental pore volume occluded for the ith point (cm³/g):

$$V_{
m inc,i} = S_{
m inc,i} R_i imes 10^{-4}$$

Cumulative pore volume occluded for the ith point (cm³/g):

$$V_{ ext{cum,i}} = V_{ ext{inc,i-1}} + V + \ldots + V_{ ext{inc,i}}$$

dV/dR pore volume for the ith point (cm³/g-Å):

$$rac{dV}{dR_{
m i}} = rac{V_{{
m inc},i}}{t_{
m i}-t_{
m i-1}}$$



QUANTITY ADSORBED FOR CHEMICAL ADSORPTION

For the ith dose:

$$N_{ ext{gas}_i} = N_{ ext{gas}_{i-1}} + V_{ ext{LOW}} imes \left[rac{P_{ ext{sys1}_i}}{T_{ ext{sys1}_i}} - rac{P_{ ext{sys2}_i}}{T_{ ext{sys2}_i}}
ight] imes rac{T_{ ext{STD}}}{P_{ ext{STD}}}$$

$$N_{ads_i} = rac{N_{gas_i} - P_{sam}}{P_{STD} imes V_{FS}}$$

$$Q_i = rac{N_{ ext{ads}_i}}{Msam}$$

where

 M_{sam} = mass of sample (g)

 N_{ads} = amount of gas adsorbed after equilibrating ith dose (cm³ STP)

 $N_{gas\ i}$ = total amount of gas dosed into sample tube after ith dose (cm³ STP)

P_{sam i} = sample pressure after equilibrating ith dose of gas onto sample (mmHg)

P_{STD} = standard pressure (760 mmHg)

P_{sys1 i} = system manifold pressure before ith dose of gas onto sample (mmHg)

P_{sys2 i} = system manifold pressure after ith dose of gas onto sample (mmHg)

 Q_i = amount of gas adsorbed per gram of sample (cm³/g STP)

T_{STD} = standard temperature

 $T_{sys1 i}$ = system manifold temperature before i^{th} dose of gas onto sample (K)

 T_{sys2} i = system manifold temperature after ith dose of gas onto sample (K)

Vfs = free space

 V_{LOW} = volume of lower manifold



QUANTITY ADSORBED FOR CHEMICAL ADSORPTION

A portion of the dosing volume may be at a slightly elevated temperature due to heating of the sample ports. The manifold volume is partitioned into a volume at the temperature of the manifold block and a volume at the average temperature of the ports.

$$egin{aligned} n_{lpha} - n_{d} - n_{f} \ & \ n_{d} = P_{1m}C(P_{1m}, T_{1m}, ar{T}_{1p}) - P_{2m}C(P_{2m}, T_{2m}, ar{T}_{2p}) \ & \ c(P, T_{m}, T_{p}) = V_{m} igg(rac{lpha}{z(P, T_{m})T_{m}} + rac{eta}{z(P, T_{p})T_{p}} igg) \end{aligned}$$

where

 α and $\beta~=~$ constants that determine the relative weights of the manifold and port tem-

peratures

n_a = quantity of gas adsorbedn_d = quantity of gas dosed

n_f = quantity of gas in the free space

P_{1m} = manifold pressure before dosing onto the sample

P_{2m} = manifold pressure after dosing

T_{1m} = manifold temperature before dosing onto the sample

T_{1p} = average of all port temperatures before dosing onto the sample

T_{2m} = manifold temperature after dosing

T_{2m} = average of all port temperatures after dosing

V_m = volume of the dosing manifold



QUANTITY ADSORBED FOR PHYSICAL ADSORPTION

FREE SPACE - MEASURED

Measured free-space volumes are calculated using the following equations:

$$V_{fw} = rac{V_{ ext{man}}}{T_{ ext{man}}}igg(rac{P_1}{P_2}-1igg)T_{ ext{STD}}$$

$$V_{fc} = V_{fw}igg(rac{P_2}{P_3}igg)$$

FREE SPACE - CALCULATED

The calculated free space is determined by subtracting the gas capacity of the volume occupied by the sample from the measured free space of the empty tube.

$$V_{fw} = V_{wb} - V_s igg(rac{T_{ ext{STD}}}{T_{ ext{amb}}}igg)$$

$$V_{
m bath} = rac{V_{fc} - V_{fw}}{1 - rac{T_{
m bath}}{T_{
m amb}}}$$

$$V_s = rac{m}{
ho}$$



COMPENDIUM OF VARIABLES

= non-ideality factor α

sample density ρ

= sample mass m

 P_i = equilibrated sample pressure

 P_1 system manifold pressure before dosing onto sample

 P_2 system manifold pressure after dosing onto sample

 P_3 sample pressure after raising dewar and equilibrating with helium

approximate room temperature (298 K) T_{amb}

T_{bath} = analysis bath temperature (K)

= system manifold temperature before dosing helium onto sample (K) T_{man}

T_{STD} = standard temperature (273.15 K)

= volume of cold free space (cm³ at standard temperature) V_{amb}

volume of cold free space at analysis bath temperature (cm³ at standard V_{bath}

temperature)

= volume of cold free space of the empty tube (cm³ at standard temperature) V_{cb}

volume of cold free space (cm³ at standard temperature) V_{fc}

volume of warm free space (cm³ at standard temperature) V_{fw}

manifold volume (cm³) V_{man}

 V_s = sample volume

= volume of warm free space of the empty tube (cm³ at standard temperature) V_{wb}



RELATIVE PRESSURE

Relative pressure is defined as:

$$P_{\mathrm{rel},i} = rac{P_i}{P_{0,i}}$$

 $P_{0,i}$ depends on the method selected in the P_0 and Temperature options

If P_0 was measured for each point, $P_{0,i}$ is the saturation pressure measured when P_i was equilibrated.

If P_0 was measured at intervals, $P_{0,i}$ is found by linear interpolation.

$$P_{0,i} = P_{0,i-1} + \left(P_{0,i+1} - P_{0,i-1}
ight) rac{t_i - t_{i-1}}{t_{i+1} - t_{i-1}}$$

The last measured P_0 is only for points taken after the last P_0 measurement.

If P_0 was calculated from measured Psat or entered temperature the calculated P_0 is used for all data points.

If P_0 was entered, the entered P_0 is used for all data points.



SATURATION PRESSURE

Saturation pressure (P_0) is selected on the P_0 and Temperature Options window. It may be entered or measured in the P_0 tube. The analyzer uses the following methods to get P_0 :

- 1. P_0 is measured in the P_0 tube for each isotherm point. This method can be used only if the instrument has a dedicated P_0 transducer.
- 2. P₀ is measured at a specified time interval using the 1000 mmHg manifold transducer.
- 3. The saturation pressure of a gas is measured in the P_0 tube for prior to data collection. The bath temperature is found by looking up the temperature for the measured saturation pressure in the Psat vs. T table. P_0 is found by looking up the calculated bath temperature in the adsorptive properties Psat vs. T table.
- 4. P₀ is found by looking up the entered bath temperature in the adsorptive properties Psat vs. T table.
- 5. P_0 is entered.



SPC REPORT VARIABLES

REGRESSION CHART VARIABLES

The line of best fit for the Regression Chart is calculated by the usual least squares method. 1 If there is only a single point or all N points have the same x-value, there can be no line of best fit in the standard form.

$$\bar{\chi} = \frac{\sum \chi_i}{N}$$

$$\bar{y} = \frac{\sum y_i}{N}$$

Slope =
$$\frac{\sum (\chi_i - \bar{\chi})(y_i - \bar{y})}{\sum (\chi_i - \bar{\chi})^2}$$

$$\text{Intercept} = \bar{y} - \text{Slope} \cdot \bar{\chi}$$

The coefficient of correlation for this line is also calculated in the usual way. 2)

$$\sigma_\chi = \sqrt{rac{\sum (\chi_4 - ar{\chi})^2}{N}}$$

$$\sigma_y = \sqrt{rac{\sum (y_i - ar{y})^2}{N}}$$

$$\mathrm{Cov}(x,y) = rac{\sum (\chi_i - ar{\chi})(y_i - ar{y})}{N}$$

$$ext{CorrelationCoeff} = rac{ ext{Cov}(x,y)}{\sigma_\chi \sigma_y}$$

¹) BASIC Scientific Subroutines Vol II, by F.R. Ruckdeschel, Copyright 1981 BYTE Publications/McGraw Hill, p. 16.

²) Mathematical Handbook for Scientists and Engineers, G.A. Korn and T.M. Korn, McGraw Hill, Sec. 18.4. (1968)



CONTROL CHART VARIABLES

$$\mathrm{Mean} = rac{\sum y_i}{N}$$

$$ext{StandardDeviation} = \sqrt{rac{\sum (y_i - ext{Mean})^2}{N-1}}$$

$$C.V. = \frac{StdDev}{Mean}$$

$$+n\sigma = \text{Mean} + n \cdot \text{StandardDeviation}$$

$$-n\sigma = \text{Mean} - n \cdot \text{StandardDeviation}$$



STOICHIOMETRY FACTOR

The calculated stoichiometry factor is a weighted average. It is dependent on both the individual stoichiometry factor and number of moles of each active metal.

$$SF_{ ext{CALC}} = rac{rac{\% ext{weight1} imes SF1}{W_{ ext{ATOMIC1}}} + rac{\% ext{weight2} imes SF2}{W_{ ext{ATOMIC2}}} + rac{\% ext{weight3} imes SF3}{W_{ ext{ATOMIC3}}} + ... + rac{\% ext{weightn} imes SFn}{W_{ ext{ATOMIC1}}}}{rac{\% ext{weight1}}{W_{ ext{ATOMIC1}}} + \% ext{weight2}W_{ ext{ATOMIC2}}} + rac{\% ext{weight3}}{W_{ ext{ATOMIC3}}} + ... + rac{\% ext{weightn} imes SFn}{W_{ ext{ATOMIC1}}}}{W_{ ext{ATOMIC1}}}$$

where

= calculated stoichiometry factor SF_{CALC} SF_1 = stoichiometry factor for first metal SF_2 = stoichiometry factor for second metal SF_3 stoichiometry factor for third metal stoichiometry factor for nth metal SF_n %weight₁ = % of sample weight for first metal %weight₂ = % of sample weight for second metal %weight₃ = % of sample weight for third metal %weight_n = % of sample weight for nth metal = atomic weight of first metal (g/mole) WATOMIC₁ WATOMIC₂ = atomic weight of second metal (g/mole) WATOMIC₃ = atomic weight of third metal (g/mole) = atomic weight of nth metal (g/mole) $WATOMIC_n$



SUMMARY REPORT

The following calculations and the results of previous calculations (as noted) are used to generate the summary report:

a. Single-point Surface Area (m²/g)

$$S_{1PT} = rac{[Q(1-P)] imes CSA(6.023 imes 10^{23})}{22414 ext{cm}^2 imes ext{STP}igg(rac{10^{18} ext{nm}^2}{ ext{m}^2}igg)}$$

where

P = pressure closest to 0.3 of the relative pressure points designated for surface area calculations

Q = quantity adsorbed corresponding to P

- b. Multi-point Surface Area. See BET Surface Area on page 2.
- c. Langmuir Surface Area. See <u>Langmuir Surface Area for Chemical Adsorption on page 41</u> and <u>Langmuir Surface Area for Physical Adsorption on page 42</u>
- d. t-Plot Micropore Surface Area. See t-Plot on page 59.
- e. t-Plot External Surface Area. See t-Plot on page 59.
- f. BJH Cumulative Adsorption
- g. BJH Cumulative Desorption
- h. Adsorption Total Pore Volume
- i. Desorption Total Pore Volume
- j. t-Plot Micropore Pore Volume. See t-Plot on page 59.
- k. Freundlich. See Freundlich Isotherm on page 29.
- I. Temkin. See Temkin Isotherm on page 60.
- m. Alpha-S. See Alpha-S Method on page 1.
- n. DFT Pore Size and DFT Surface Energy. See <u>DFT (Density Functional Theory) on page 14</u>.
- o. Nanoparticle Size

$$d = rac{6 imes 10^4}{A
ho}$$

where

 ρ = sample density

A = BET surface area

d = side length (for cubic particles or diameter (for spherical particles)

- p. Dubinin-Astakhov Micropore Surface Area. See **Dubinin-Astakhov on page 19**.
- q. Dubinin-Astakhov Micropore Volume. See Dubinin-Astakhov on page 19.



- r. Dubinin-Radushkevich Micropore Surface Area. See <u>Dubinin-Radushkevich on page 23</u>.
- s. Dubinin-Radushkevich Monolayer Capacity. See Dubinin-Radushkevich on page 23.
- t. MP-Method Cumulative Surface Area of Pores

$$S_{
m total} = S_{
m cum,i}$$

See <u>MP-Method on page 46</u> for the last collected data point used in the MP-method Calculations, and the range of hydraulic pore radii over which the cumulative surface area was computed.

u. MP-Method Cumulative Pore Volume of Pores

$$V_{
m total} = V_{
m cum,i}$$

See <u>MP-Method on page 46</u> for the last collected data point used in the MP-method calculations, and the range of hydraulic pore radii over which the cumulative pore volume was computed.

v. Average Pore Hydraulic Radius (Å)

$$ar{r} = rac{V_{
m total}}{S_{
m total}} imes 10^4$$

w. Horvath-Kawazoe. See Horvath-Kawazoe on page 31.



T-PLOT

A least-squares analysis fit is performed on the $(t_i, N_{\mathsf{ads},i})$ data pairs where t_i is the independent variable and $N_{\mathsf{ads},i}$ is the dependent variable. Only the values of t_i between t_{min} and t_{max} , the minimum and maximum thickness, are used. The following are calculated:

- a. Slope (S cm³/g-Å STP)
- b. Y-intercept (Y_{int} cm³/g STP)
- c. Error of the slope (S_{err} cm³/g-Å STP)
- d. Error of the Y-intercept (YI_{err} cm³/g STP)
- e. Correlation coefficient

Using the results of the above calculations, the following can be calculated:

External Surface Area (m²/g):

$$\frac{\mathit{SV}_{\mathrm{mol}}}{\mathit{F} \times 22414\mathrm{cm}^3\mathit{STP}} \times 10^4$$

where

 10^4 = unit conversions

F = surface area correction factor, user-entered on the *t-Plot Report Options* window

 V_{mol} = liquid molar volume, from the fluid property information

Micropore Surface Area (m²/g):

$$SA_{\mu
ho} = SA_{
m total} + SA_{
m ext}$$

where SA_{total} is the BET surface area if the user enabled the BET report exclusively, or Langmuir surface area if the user enabled the Langmuir report exclusively. If neither report has been selected, SA_{total} is the BET surface area value calculated using a set of default parameters.

Micropore Volume (cm³ liquid/g):

$$rac{Y_{
m int}V_{
m mol}}{22414{
m cm}^3{
m STP}}$$



TEMKIN ISOTHERM

The Temkin isotherm has the form:

$$rac{Q}{Q_{m}}=rac{RT}{q_{0}lpha} ext{ln}(A_{0}P)$$

where

 A_0 = adjustable constant

a = adjustable constant

P = equilibrium pressure measured by gauge at temp T_{amb}

 q_0 = the differential heat of adsorption at zero surface coverage

Q = quantity of gas adsorbed

 $Q_{\rm m}$ = quantity of gas in a monolayer

R = molar gas constant $8.31441 \times 10^{-3} \frac{kJ}{\text{mol}K}$

T = bath temperature

In terms of quantity adsorbed

$$Q=rac{RTQ_m}{q_0lpha}[\ln A_0+\ln P]$$

Thus, the plot of the natural log of absolute pressure vs. quantity adsorbed yields a straight line with slope $\frac{RTQ_{\rm m}}{q_0}$ and intercept $\ln A_0 \frac{RTQ_{\rm m}}{q_0 \alpha}$



THERMAL TRANSPIRATION CORRECTION

During data reduction, thermal transpiration correction is applied to the data if the user selected *Apply thermal transpiration correction* from the *Report Options* window. Starting with the first collected pressure, the following calculations are performed until the pressure ratio (PC/P) is greater than or equal to 0.99.

$$Y = \left(rac{P imes ext{SD} imes ext{MD}^2}{2.33 imes T}
ight)\! 10^3$$

$$\mu = \frac{(1+G)Y}{(1+H)Y}$$

$$F=rac{1}{lpha Y^2+eta Y+\mu}$$

$$P = \left(1 - Figg(1 - \sqrt{rac{T_{
m bath}}{T_{
m amb}}}igg)
ight)$$

where

 α = Weber's coefficient, 0.033

 β = Weber's coefficient, 0.245

 F, Y, μ = intermediate values for subsequent calculations

G = Weber's coefficient, 2.5 H = Weber's coefficient, 2

MD = thermal transpiration hard sphere diameter of gas (Å), from the Adsorptive Prop-

erties window

P = equilibrated collected pressure measured by gauge at temp T_{amb}

SD = inside diameter of sample tube (mm) from the Report Options window

T = average temperature $\frac{T_{\text{bath}} + T}{2}$

 T_{amb} = room temperature (298 K)

 T_{bath} = analysis bath temperature (K), from the P_0 and Temperature Options window



THICKNESS CURVE

For each point designated, the following parameters are used in thickness curve calculations:

C₁ = parameter #1

C₂ = parameter #2

 C_3 = parameter #3

 $P_{rel, i}$ = relative pressure for the ith point (mmHg)

t_i = thickness for ith point

REFERENCE

Interpolated from table.

KRUK-JARONIEC-SAYARI

$$t_{
m i} = \left(rac{C_2}{C_2 - \log(P_{{
m rel},i})}
ight)^{c^3}$$

HALSEY

$$t_{ ext{i}} = C_1 igg[rac{C_2}{\ln(P_{ ext{rel,i}})}igg]^{C_3}$$
 Halsey 1)

HARKINS AND JURA

$$t_{
m i} = \left\lceil rac{C_1}{C_2 - \log(P_{
m rel,i})}
ight
ceil^{C_3}$$
 Harkins and Jura 2)

¹⁾ Halsey, G., J. Chem. Phys. 16, 931-937 (1948).

²⁾ Harkins, W.D. and Jura, G., J. *Chem. Phys.* 11, 431 (1943).



BROEKOFF-DE BOER

$$\log(P_{\mathrm{rel,i}}) = rac{C_1}{t^2,\mathrm{i}} + C_2 \exp(c_3 t_\mathrm{i})$$

CARBON BLACK STSA

$$t_{
m i} = C_1 (P_{
m rel,i})^2 + C_2 (P_{
m rel,i}) + C_3$$



Blank Page